# BAMBANKER™

## DMSO Free

DMSO-free and Serum-free  
cell freezing medium

<For Research Use Only>

**BAMBANKER™** DMSO Free is a DMSO-free and serum-free medium and preservation of valuable culture cells such as tumor cells.

### Features

- <1> Ready-to-use medium for preservation of cells, <2> Usable without diluting, <3> No program freezer is required, <4> DMSO-free and Serum-free

### Operating Procedure

- 1) Centrifuge the cells in logarithmic growth phase, and collect  $5 \times 10^5 \sim 1 \times 10^7$  cells.
- 2) Suspend the collected cells in 1 mL of this medium, and place the cells in cryotubes for freezing and preservation. Then, freeze and preserve the cells at  $-80^\circ\text{C}$  without preliminary freezing. The cells frozen at  $-80^\circ\text{C}$  can be preserved subsequently in liquid nitrogen.
- 3) Thawing must be carried out quickly in a constant-temperature chamber or bath.
  - ★ The cells must be frozen in logarithmic growth phase.

### Sterility test

- Endotoxin: Colorimetric method  
Mycoplasma: Culture method  
Fungi and bacteria: Agar streak test

### Applications <Cells preserved well by cryopreservation test>:

- Jurkat (Human leukemia T cell line)  
HEK293 (Human embryonic kidney cell line)  
HepG2 (Human liver cancer cell line)

### Precautions

- ◆ Use only for research purposes, not for human use.
- ◆ Prior to using this medium, perform the confirmation test on the cells under study.
- ◆ The manufacturer shall not be responsible for any accident or damage caused by the use of this product.
- ◆ If you have any questions on the use of this medium, please contact the local distributor.

Genetics Catalog No.	Description	Package Size	Storage	Expiration period
BBF01	DMSO-free Serum-free Cell Freezing Medium	20mL	Keep at 2~10°C	Two years from manufacturing

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